

IPTG (Dioxane-Free)

Shipping: On Dry/Blue Ice

Catalog numbers

Batch No.: See vial

BIO-37036 : 5 g



A Meridian Life Science® Company

Store at -20°C

Storage and stability:

The IPTG is shipped on dry/blue ice. On arrival store at -20 °C for optimum stability. Repeated freeze/thaw cycles should be avoided.

Expiry

When stored under the recommended conditions and handled correctly, full activity of the kit is retained until the expiry date on the outer box label.

Safety precautions: Harmful if swallowed. Irritating to eyes and skin. Please refer to the material safety data sheet for further information and safety handling practice.



Xi: IRRITANT

Quality control specifications:

IPTG is extensively tested for activity and absence of contamination prior to release.

Notes:

Research use only.

Description

Isopropyl-β-D-thiogalactopyranoside (IPTG) is a chemical analogue of galactose, which cannot be hydrolysed by the enzyme β-Galactosidase. Hence, it induces the *E. coli lac* operon activity by binding and inhibiting the lac repressor without being degraded. Genes controlled by the *lac* or *tac* promoter/operator sequences are expressed to high levels in the presence of IPTG.

It is recommended for blue/white screening by *lacZ* α-complementation with appropriate vectors and host strains, such as alpha-select, CH3-Blue and *BIOBlue* cells, and for protein expression with BL21 (DE3) cells and other lac promoter-derived expression systems.

Features

- Induces *E. coli lac* operon activity
- >99.6% purity by HPLC
- Available as powder and stabilized stock solution

Applications

- Blue/white color screening
- Induction of *lac* operon for protein expression
- Genes controlled by the *lac* or *tac* promoter/operator sequences are expressed to high levels in the presence of IPTG

Characteristics

Empirical formula	C ₉ H ₁₈ O ₅ S
Molecular weight	238.3
Form	1M Solution in water
Melting point	112-113 °C
Purity	≥99% by HPLC

Recommendations for use

- Dissolve to desired concentration in Water, 18.2MΩ (DNase/RNase-free).
- Filter sterilize through a 0.22 micron disposable filter.
- Store solution at -20 °C.
- For blue/white screening, 0.1mM final IPTG concentration in LB (Luria Broth) media is recommended

Associated Products

Product Name	Pack Size	Catalog No.
X-GAL Powder	1 g	BIO-37035
Quick-Stick Ligase	50 reactions	BIO-27027
SureClean Plus	5 ml	BIO-37047
α-Select Gold Efficiency Competent Cells	1 ml	BIO-85027
CH3-Blue 10 ⁹ Chemically Competent Cells	1 ml	BIO-85040
BIOBlue 10 ⁹ Chemically Competent Cells	1 ml	BIO-85037

Product Citations:

1. Ali, F.R., *et al. Meth. Mol. Biol.* **628**, 195-214 (2010).
2. Prabhakar, V., *et al. FEBS Lett.* **583(6)**, 983-991 (2009).
3. Chan, C-H., *et al. Conserv. Gen.* **9(4)**, 1067-1070 (2008).
4. Hamblin, K., *et al. Mol. Microbiol.* **68(6)**, 1395-1405 (2008).
5. Maruta, F., *et al. J. Drug Targeting* **15(4)**, 311-319 (2007).
6. Ross, P.J., *et al. Infect. Immun.* **72(3)**, 1568-1579 (2004).
7. Meng, G. & Fütterer, K. *Nat. Struct. Biol.* **10**, 935-941 (2003).

Bioline Reagents Ltd
UNITED KINGDOM
Tel: +44 (0)20 8830 5300
Fax: +44 (0)20 8452 2822

Bioline USA Inc.
USA
Tel: +1 508 880 8990
Fax: +1 508 880 8993

Bioline GmbH
GERMANY
Tel: +49 (0)337 168 1229
Fax: +49 (0)3371 68 1244

Bioline (Aust) Pty. Ltd
AUSTRALIA
Tel: +61 (0)2 9209 4180
Fax: +61 (0)2 9209 4763

Bioline France
FRANCE
Tel: +33 (0)1 42 56 04 40
Fax: +33 (0)9 70 06 62 10

Meridian Bioscience Asia Pte Ltd
SINGAPORE
Tel: +65 6774 7196
Fax: +65 6774 6441