

Isolation of fecal DNA



Vortex Bashing Beads Lysis
Tube with fecal or soil sample



10,000 x g, 1 min



Filter lysate with Spin II A
Filter



8000 x g, 1 min



Bind, Wash, Elute DNA with
Spin Column



10,000 x g, 1 min



Filter DNA with Spin II B Filter



8000 x g, 3 min



Transfer eluted DNA



16,000 x g, 3 min



Isolated DNA

Technical Support

The troubleshooting guide is provided in the online protocol.
For technical support please email us at tech@bioline.com
or visit www.bioline.com